


Dear Phywe customer,

We congratulate you on your selection of our software *measure for Cobra4*. The combination of the Cobra4 interface system with software *measure* opens up a magnitude of possibilities for data recording and analysis.

Installation of the software


Insert the DVD into the DVD drive. As a rule, installation starts automatically. Select the menu prompt "Install or update *measure*" , to install *measure for Cobra4*. At the end of the installation process the Activation Assistant is opened. Choose "activate" and enter the enabling code that is printed on the inside of the DVD jacket. With this, the Cobra4 functionality of the software *measure* is enabled. The Activation Assistant is always available in *measure* in menu "Options" to enable further devices, if necessary.


Should a version of *measure* be already installed on your computer, this will be automatically recognized and an update of the available components will be carried out.


When the latest version of *measure* is running on your computer, all previous versions can be uninstalled.

Putting the system into operation


Details on the handling of Cobra4 instruments (USB-Link and Wireless-Link) are to be found in the operating instructions supplied with them.

To open *measure for Cobra4*, start the software *measure*  after you have installed the full version and entered the enabling code for Cobra4 as described above.

There are two ways to open *measure for Cobra4*: (1) In *measure* either via the menu button  resp. via the selection of "Cobra4" in menu „Gauge“, or (2) automatically by connecting one or several Cobra4 instruments. Connect your Cobra4 USB-Link or Wireless-Link with a Sensor-Unit plugged in as described in the operating instructions. When the drivers are installed, the display of the Cobra4 USB-Link resp. Wireless-Link lights up and the programme Navigator shows the corresponding Cobra4 instrument with connected Sensor-Unit.

You can now measure! As suggestions we have included descriptions of experiments in the programme as pdf documents that can be called up with the  button or in the menu "Experiment → Open experiment" (for this, Adobe Reader 8, also included on the installation DVD, must be installed). The necessary settings for carrying out measurements and the measurement examples are loaded automatically as well.

A complete set of all presently available experimental protocols is included in the manual "Computer-assisted experiments with Cobra4". This manual can be ordered from PHYWE using item no. 01330.02. How to carry out a measurement is described in detail in the appropriate experimental protocol together with context-related help topics that provide further information on the programme functions.

You can also call up help texts on various themes, such as "First steps with Cobra4" via button  or the "Help" menu.

If you have any questions or need further assistance, please do not hesitate to contact us.

We hope you will enjoy using Cobra4!