

Image construction at the camber mirror



Physics

Light & Optics

Reflection & refraction of light



Difficulty level

easy



Group size

2



Preparation time

10 minutes




Execution time

10 minutes

PHYWE
excellence in science

Teacher information

Application

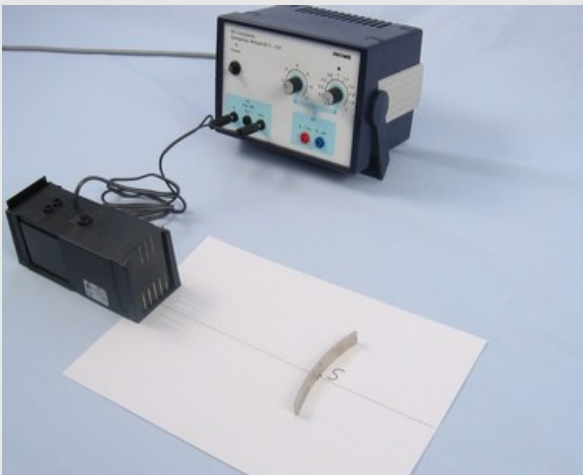PHYWE
excellence in science

Reflection at the camber mirror

A convex mirror is a mirror that is curved outwards. Usually parts of spheres are used for this purpose, onto which the light falls from the outside. Especially in road traffic, such mirrors are used to view larger areas or unclear places.

But also at Christmas time we encounter Wöb Spiegel in the form of shiny Christmas tree baubles. If you look at the mirror image, you can see a reduced, upright image.

Other teacher information (1/4)

PHYWE
excellence in science

Previous



The students should have previously learned the basics of the rectilinear propagation of light and the application of the law of reflection. They should also be able to draw individual ray paths and construct pixels.

Principle



Convex mirrors have a characteristic beam path. The reflected light rays appear to come from a point behind the mirror. Light beams incident on the convex mirror parallel to the optical axis run outwards after reflection, they diverge. Light rays incident along the optical axis are reflected within themselves.

Other teacher information (2/4)

PHYWE
excellence in science

Learning



With this experiment, the students are to become acquainted with a procedure that makes it possible to construct the image on the convex mirror for a given object. Selected light beams and their characteristic course are used for this purpose.

Tasks



The aim of this experiment is the experimental determination of the intersection point of selected light beams incident on a convex mirror and the resulting possibility of image construction. In a preliminary experiment, the properties of the mirror image are first determined, thus making it possible to compare the two experimental results later.

Other teacher information (3/4)

PHYWE
excellence in science

Additional information

The experiment is very demanding in terms of the assumed experimental skills of the students and the knowledge that is gained. No real images are obtained. The intersection point of the backward extended reflected light beams lies behind the convex mirror.

This gives the student the opportunity to explain the nature of virtual (apparent) images and to deepen his knowledge of virtual images on the plane mirror.

Other teacher information (4/4)

PHYWE
excellence in science

Notes on structure and implementation

It must be ensured that the adjustment of the mirror (centre of the outward curved surface is at point S of the optical axis) and the light box (incidence of the narrow beam of light along the optical axis to check the correct setting) is carried out very carefully by the student in order to arrive at a clear and convincing result.

To facilitate the parallel displacement of the light box, a thin pencil line parallel to the optical axis should be drawn through the tip of the object arrow to the camber mirror before the experiment.

Safety instructions

PHYWE
excellence in science

The general instructions for safe experimentation in science lessons apply to this experiment.

PHYWE
excellence in science

Student Information

Motivation

PHYWE
excellence in science


We encounter mirrors every day in a wide variety of designs. A special type of mirror is the so-called convex mirror. This is a mirror that is curved outwards. Often, partial surfaces of spheres are used for this purpose, onto which the light shines from the outside.

A typical example of a convex mirror is a Christmas tree ball, as shown in the picture on the right. Have you ever looked at your reflection in a sphere, do you notice anything special?

Christmas tree balls

Task

PHYWE
excellence in science

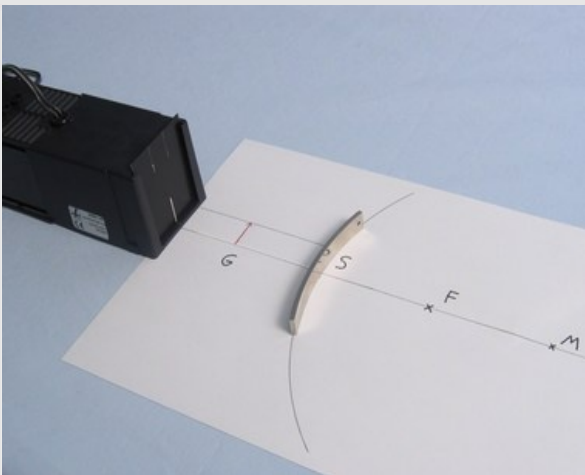

Test setup

Why do you see yourself diminished in a shiny Christmas bauble?

1. investigate the properties of the mirror on a convex mirror.

Investigate the formation of images on a convex mirror using selected light beams.

Material


Position	Material	Item No.	Quantity
1	Light box, halogen 12V/20 W	09801-00	1
2	Mirror, concave-convex	09812-00	1
3	PHYWE Power supply, 230 V, DC: 0...12 V, 2 A / AC: 6 V, 12 V, 5 A	13506-93	1

Additional material

PHYWE
excellence in science

Position	Material	Quantity
1	Ruler (approx. 30cm)	1
2	White paper (DIN A4)	1
3	Circle	1

Structure

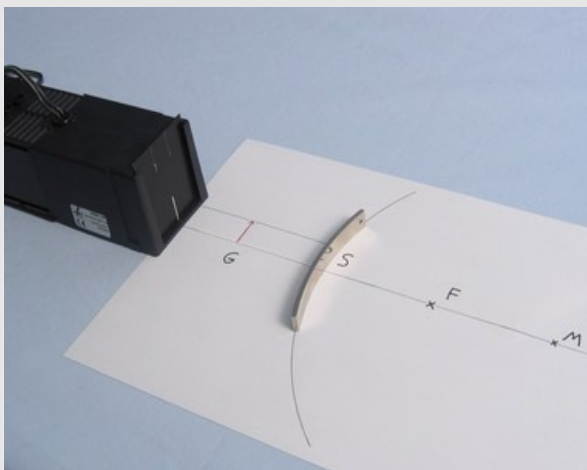PHYWE
excellence in science

Test setup

Attention


Make sure that the convex mirror is always aligned with the center of the outer camber on the point S and does not change its position when the light box is moved.

1. properties of the mirror image

- Look into the convex mirror held vertically in front of your face. What characteristics of your reflection do you notice? Write down your observations in the protocol.

Procedure (1/5)

PHYWE
excellence in science

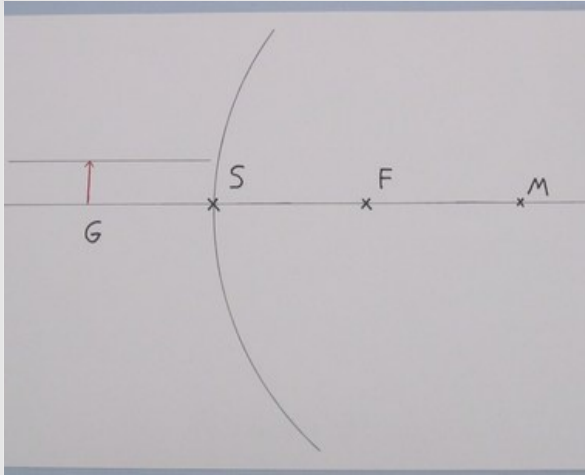

Moving the light box



2. image formation at the camber mirror

- Prepare a sheet of paper as shown on the left
- The routes \overline{FS} and \overline{MS} are 7.2 cm each; the circle around M has the radius \overline{MS} .
- Draw at a distance of 6 cm from the point S (opposite of F and M) draw a vertical arrow, 2 cm long, on the optical axis and denote it by G (Subject).
- Draw a thin auxiliary line parallel to the optical axis, passing through the tip of the arrow.

Procedure (2/5)


PHYWE
excellence in science

- Insert the slit diaphragm into the light box on the lens side. Place the light box and the convex mirror on your sheet of paper.
- Connect the light box to the power supply (12 V ~).
- Check the adjustment of the camber mirror by letting the narrow beam of light fall along the optical axis.

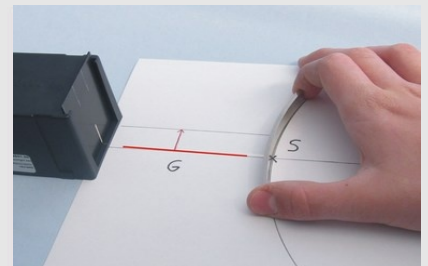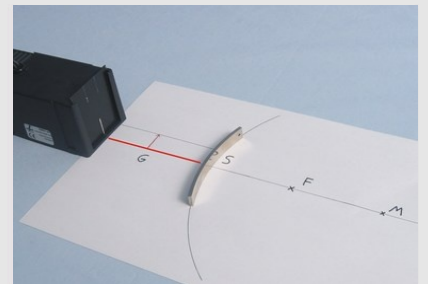

Implementation (3/5)

PHYWE
excellence in science

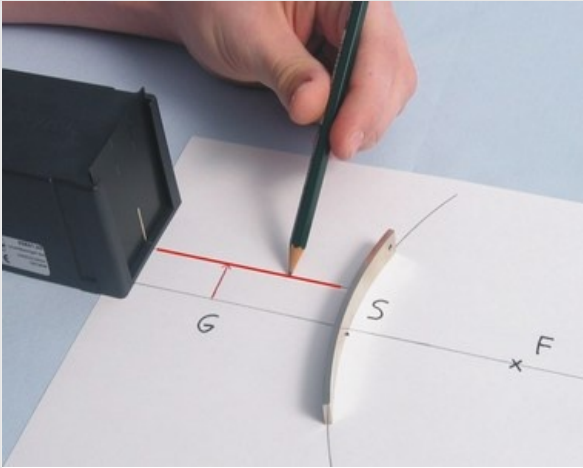


Moving the light box

- Now move the light box until the narrow light beam runs parallel to the optical axis along the auxiliary line and through the tip of the arrow (imaginary object).
- Observe the light beam reflected by the convex mirror and mark the course of the incident and reflected light beam with two crosses each.

Procedure (4/5)

PHYWE
excellence in science

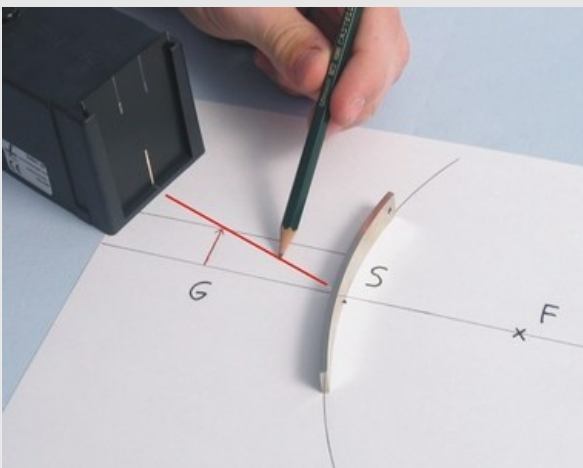

Turning the light box

- Turn the light box until the light beam is exactly in the direction of the dot S but still through the tip of the object arrow G is running.
- Again, observe the light beam reflected by the convex mirror and mark the course of the incident and reflected light beam twice with crosses (use a different color or marker).

Procedure (5/5)

PHYWE
excellence in science



Picture of the power supply unit

- Turn off the power supply and remove the light box and the camber mirror from the sheet of paper.
- Connect the matching crosses so that the course of the light beams before and after the reflection at the convex mirror becomes clear.
- Extend the reflected light rays beyond the mirror with broken lines. What can you determine?
- Write down your observations.

PHYWE
excellence in science



Protocol

Task 1

Consider the following statement:

If the object is at a distance of 6 cm from the convex mirror, the image is at a shorter distance but on the other side.

The image is reduced and upright (it is a virtual/apparent image).

 True Incorrect Check

Task 2

Why do you see yourself diminished in a shiny Christmas bauble?

A [] Christmas tree ball is a small []. Thus, the [] and the laws about the course of selected [] also apply to it.

As a result, a [], upright, (apparent) [] results from the [].

shiny

observer

image

curved mirror

diminished

light beams

law of reflection

 Check

Task 3

PHYWE
 excellence in science

Draw a straight line from the tip of the object arrow G to the point M . What can you tell?

How would a beam of light incident on the convex mirror along this straight line be reflected?

The also passes through the of the other two selected light rays and thus through the of the image arrow on its way to the lying behind the concave mirror. A light beam incident on the behind the concave mirror is reflected into itself (angle of incidence = =).

Task 4

PHYWE
 excellence in science


What is special about the images created at the camber mirror?

The images on the convex mirror are behind the mirror. Therefore, as with the plane mirror, they are apparent images.

The images on the camber mirror are in front of the mirror. They are therefore real images

Slide	Score/Total
Slide 20: Properties of the image	0/1
Slide 21: Christmas ball as a camber mirror	0/7
Slide 22: Interpretation of the beam path	0/7
Slide 23: Apparent images at the camber mirror	0/1

Total  0/16

 Solutions

 Repeat